

MIKRO 220 | R

MICRO / GENERAL PURPOSE / LIPEMIC

High RCF, high-throughput and highly versatile for time saving performance. The MIKRO 220 | R features a high capacity 48-place angle rotor and 60-place drum rotor for maximum productivity. Higher RCF values reduce production time. Additional rotors include a 24-place swing-out rotor for microliter tubes, 6-place PCR strip rotor and rotors for conical tubes, blood tubes, pediatric tubes, urine tubes and hematocrit capillaries.

Robust and quick temperature control saving time and energy. Available with optional temperature control (-20 to +40 °C), the MIKRO 200 R cools down to +4°C in just 10 to 15 minutes using the Fast-Cool function.

Programmable for easy recall of application parameters limiting user error. The flexible control panel further enables simplified and secure centrifugation by providing up to 10 user-defined programs for easy recall of adjustable application parameters.

Low cost alternative for Lipemic samples. The MIKRO 220 eliminates the need for a limited use, high-cost, ultra-centrifuge and offers a solution for clearing Lipemic samples at a fraction of the cost. Process your samples quickly, in-house, and do away with send-outs.



MODEL	MIKRO 220	MIKRO 220 R
Max. Capacity	48/60 x 1.5 - 2.0 ml	48/60 x 1.5 - 2.0 ml
Max. RPM / RCF	18,000 / 31,514	18,000 / 31,514
Max. Blood Tubes*	12 (up to 10 ml)	12 (up to 10 ml)
Max. Conical Tubes*	6/3 (15/50 ml)	6/3 (15/50 ml)
Temperature Control	Ambient	-20 °C to +40 °C
Dims (H x W x D)	12 x 13 x 16.5 in.	12 x 13 x 25.5 in.
Warranty [†]	5 yr.	5 yr.

*Please inquire about the fit of your specific tube dimensions. † Warranty valid in U.S. and Canada only, see warranty statement for complete details

Learn more: hettweb.com/mikro220



MIKRO 220 | R CENTRIFUGE PACKAGES

MIKRO 220 MICROLITER TUBE PACKAGE - 24 - BIO-CONTAINMENT



Tube Qty.	Volume (mL)	Size (mm)	RPM	RCF	
24	1.5 - 2.0		≤ 18,000	≤ 31,514	
Includes: (1) Centrifuge (1) Fixed-angle Rotor (24) Adapters 1.5 ml (1) Rig-containment Lid					

PKG. NO.	PKG. NO. (TEMP. CONTROL)
220MICR024	220RMICR024

MIKRO 220 MICROLITER TUBE PACKAGE - 30 - BIO-CONTAINMENT



Tube Qty.	Volume (mL)	Size (mm)	RPM	RCF	
30	1.5 - 2.0		≤ 14,000	≤ 21,255	
Includes: (1) Centrifuge,	Includes: (1) Centrifuge, (1) Fixed-angle Rotor, (30) Adapters 1.5 mL, (1) Bio-containment Lid				

PKG. NO.	PKG. NO. (TEMP. CONTROL)
220MICR030	220RMICR030

MIKRO 220 MICROLITER TUBE PACKAGE - 48 - BIO-CONTAINMENT



Tube Qty.	Volume (mL)	Size (mm)	RPM	RCF
48	1.5 - 2.0		≤ 14,000	≤ 18,845
Includes: (1) Centrifuge,	(1) Fixed-angle Rotor, (48)	Adapters 1.5 mL, (1) Bio-c	ontainment Lid	

PKG. NO.	PKG. NO. (TEMP. CONTROL)
220MICR048	220RMICR048

MIKRO 220 MICROLITER TUBE PACKAGE - 90 - BIO-CONTAINMENT



Tube Qty.	Volume (mL)	Size (mm)	RPM	RCF
24	1.5 - 2.0		≤ 13,000	≤ 18,516
Includes: (1) Centrifuge,	(1) Horizontal Rotor, (24) A	dapters 1.5 mL, (1) Bio-cor	ntainment Lid	

PKG. NO.	PKG. NO. (TEMP. CONTROL)	
220MICR090	220RMICR090	

MIKRO 220 MICROLITER TUBE PACKAGE - STRIP



Strip Qty.	Volume (mL)	Size (mm)	RPM	RCF
6	PCR Strip: (8) 0.2		≤ 14,000	≤ 18,845
Includes: (1) Centrifuge,	(1) Fixed-angle Rotor			

PKG. NO.	PKG. NO. (TEMP. CONTROL)
220MICROSTRIP	220RMICROSTRIP

MIKRO 220 CRYO TUBE PACKAGE



Tube Qty.	Volume (mL)	Size (mm)	RPM	RCF
20	1.8		≤ 14,000	≤ 18,407
Includes: (1) Centrifuge,	(1) Fixed-angle Rotor, (1) B	lio-containment Lid (optiona	al)	

	PKG. NO.	PKG. NO. (TEMP. CONTROL)	
	220CRY0	220RCRY0	
	For bio-containment add "-BC" to package no. above		

MIKRO 220 R CELL CULTURE PACKAGE



Tube Qty.	Volume (mL)	Size (mm)	RPM	RCF
6	15	17 x 120	≤ 6,000	≤ 3,824
3	50	29 x 115	≤ 6,000	≤ 3,824
Includes: (1) Centrifuge, (1) Fixed-angle Rotor, (6) Adapters 15mL, (3) Adapters 50 mL				

PKG. NO). (TEMP. CONTROL)
220RCE	LLCULTURE

MIKRO 220 CUSTOM PACKAGE



Supports over 82 application configurations, please inquire for more information

Tube types include: Microliter, Cryo, PCR strips, Pediatric, Blood, Standard and Conical up to 50 mL, Hematology

Learn more:	hettweb.com/	mikro220/
-------------	--------------	-----------

Hettich | North America

100 Cummings Center, Ste. 136L Beverly, MA 01915, U.S.A. Toll free: 1 (866) 370-4388 info@hettweb.com

hettweb.com

© 2019 Hettich Instruments, LP. All rights reserved. Lit. No. PS-MIKRO-220-EN-01-2019.

Hettich products are designed to help you achieve optimal results for your application and are built to perform to the specifications outlined in the operators manual. For application specific information and settings, please refer to your organizations standard operating procedure. As always, our Hettich representatives are here to help determine which Hettich products and/or accessories best fit your laboratory requirements.